

QUANTOFIX® Sulfate

en

Pack content:

1 aluminum container with 100 test strips

Measuring range:

200–1600 mg/L SO_4^{2-}

Color gradation:

< 200 · > 400 · > 800 · > 1200 · > 1600 mg/L SO_4^{2-}

General indications:

Remove only as many test strips as are required. Close the container immediately after removing a strip. Do not touch the test fields.

Additional indications:

If the sample solution has a pH value of 4–8, the reaction will take place without interferences. Strongly acid solutions have to be buffered with sodium acetate, and alkaline solutions with tartaric acid.

Instructions for use:

1. Dip the test strip with all test fields into the sample for 1 s.
2. Shake off excess liquid.
3. Wait 120 s.
4. Compare test field with the color scale.

Interferences:

The following ions interfere with the reading only above the indicated concentrations:

- > 1000 mg/L BrO_3^- , ClO_3^- , CrO_4^{2-} , SO_3^{2-} , $\text{S}_2\text{O}_5^{2-}$, SeO_3^{2-} ,
- > 500 mg/L $\text{S}_2\text{O}_4^{2-}$,
- > 100 mg/L CN^- , S_2^- .





Disposal:







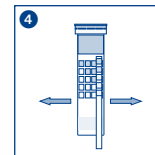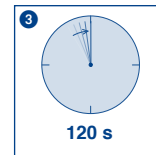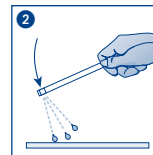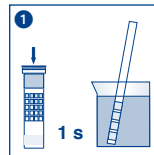


Used test strips can be placed in the normal household waste

Storage:

Avoid exposing the strips to sunlight and moisture. Keep container cool and dry (storage temperature not above +30 °C).

If correctly stored, the test strips may be used until the use-by-date printed on the packaging.



 REF	Artikelnummer / Item number / Référence produit / Referencia / Artikelnummer / Numero di catalogo / Numer artykułu	 LOT	Chargencode / Batch identification / Número de lot / Código de lote / Produktionummer / Codice del lotto / Numer partii
	Achtung! / Warning! / Attention! / ¡Atención! / Let op! / Attenzione! / Uwaga!		Verwendbar bis / Use by / À utiliser avant / Fecha de caducidad / Te gebruiken tot / Utilizzare entro (anno / mese) / Przydatność do użycia
	Temperaturbegrenzung / Permitted storage temperature range / Limites de température / Limites de temperatura / Temperatuurbegrenzing / Limite de temperatura / Ograniczenie temperatury		Trocken aufbewahren / Store in a dry place / Conserver au sec / Mantener seco / Op een droge plaats bewaren / Mantenerse asciutto / Przechowywać w suchym miejscu
	Gebrauchsanweisung beachten / Please read instructions for use / Respecter les instructions d'utilisation / Observarse las instrucciones de uso / Lees de bijsluiter / Consultare le istruzioni per l'uso / Przestrzegać instrukcję użycia		Packung geschlossen halten / Keep container closed / Refermer la boîte / Mantenga el envase cerrado / Verpak- king gesloten houden / Conservare la confezione chiusa / Przechowywać pojemnik szczelnie zamknięty
	Ausreichend für <n> Prüfungen / Contains sufficient for <n> tests / Contenu suffisant pour <n> tests / Contenido suficiente para <n> tests / Voldoende voor <n> tests / Contenuito sufficiente per <n> test / Wystarczająco dla kontroli <n>		