# visocolor<sup>®</sup> alpha

## Nitrat

### Nitrate / Nitrati / Nitrato / Nitraat

ACHTUNG  
WARNING  
ATTENTION

REF 935 065

2–50 mg/L NO<sub>3</sub><sup>-</sup>

100 Tests

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#### Test kit for performing colorimetric tests on nitrate

en

**Hazard warning:** NO<sub>3</sub>-1 contains acetic acid 0–25%, NO<sub>3</sub>-2 contains sulfanilic acid 80–100%.

H317 May cause an allergic skin reaction.

**Instructions for use:** also refer to the pictogram on the inside

1. Rinse the test vessel several times with the water sample and fill to the ring mark (5 mL).
2. Add 5 drops of NO<sub>3</sub>-1 and mix by swirling.
3. Add 1 level measuring spoon of NO<sub>3</sub>-2 and swirl for 30 s.
4. After 5 min place the measuring vessel on the color chart and assign the value by comparison of the color. Mid-values can be estimated.

This method can also be used for analyzing sea water.

**Disposing of the sample:** The used analysis specimens can be flushed down the drain with tap water and channelled off to the local sewage treatment works.

**Interferences:** Nitrate concentrations > 50 mg/L NO<sub>3</sub><sup>-</sup> lead to an instable color reaction, the reaction color gets bleached. Dilute sample until the concentration is within the measuring range. Nitrite interferes. Depending on their concentration, oxidizing substances, such as Cl<sub>2</sub>, Cr(VI) or Fe(III), may reduce the measurement reading or suppress the reaction totally.

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