

Code 3031

High Range
Free Chlorine


50 Test Strips



Insta-Test[®] ANALYTIC

High Range Free Chlorine					
■	■	■	■	■	■
0	50	100	250	500	800

Using a cup-size sample, immerse strip for **2 seconds**. Remove strip with **pad face up**. **DO NOT SHAKE OFF EXCESS WATER**. With pad still facing up. Immediately compare to color chart. Discard strip after 30 seconds.

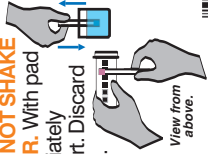

- Keep wet fingers out of the bottle.
- Cap bottle tightly.
- Store in a cool, dry place.

LaMotte Company
Chesterown, MD 21620
<http://www.lamotte.com>

